Colisure^{*}

06-14048-06

IDEXX

For Technical Support, please call:

North/South America: 1 207 556 4496/1 800 321 0207

Europe: 00800 4339 9111

UK: +44 (0) 1638 676800

China: +86 21 61279528

Japan: +03 5301 6800

Australia: 1300 443 399

IDEXX

IDEXX Laboratories, Inc., One IDEXX Drive, Westbrook, Maine 04092 USA

idexx.com/water

Colisure* Test Kit

Introduction

Colisure* simultaneously detects total coliforms and *E. coli* in water. It is based on IDEXX's proprietary Defined Substrate Technology*. When total coliforms metabolize Colisure's DST* nutrient-indicator, CPRG, the sample turns from yellow to red/magenta. When *E. coli* metabolize Colisure's DST* nutrient-indicator, MUG, the sample also fluoresces. Colisure can simultaneously detect these bacteria at 1 cfu/100 mL within 24 hours even with as many as 2 million heterotrophic bacteria per 100 mL present.

Storage

Store at 2–25°C away from light.

Presence/Absence (P/A) Procedure

1. Allow sample to equilibrate to room temperature.
2. Add contents of one pack to a 100 mL sample in a sterile, transparent, nonfluorescing vessel.
3. Cap vessel and shake.
4. Incubate at 35±0.5°C for 24 hours.
5. Read results according to Result Interpretation Table below.

Quanti-Tray* Enumeration Procedure

1. Add contents of one pack to a 100 mL water sample in a sterile vessel.
2. Add 5 drops of IDEXX Antifoam Solution (Catalog # WAFDB).
3. Cap vessel and shake thoroughly. Allow sample to stand for at least one minute and shake thoroughly again to make sure there are no large media particles remaining.
4. Pour sample/reagent mixture into a Quanti-Tray* or Quanti-Tray*/2000 and seal in an IDEXX Quanti-Tray* Sealer.
5. Place the sealed tray in a 35±0.5°C incubator for 24 hours.
6. Read results according to the Result Interpretation table below. Count the number of positive wells and refer to the MPN table provided with the trays to obtain a Most Probable Number.

Result Interpretation

Appearance	Result
Yellow/gold	Negative for total coliforms and <i>E. coli</i>
Red or magenta	Positive for total coliforms
Red/magenta and fluorescence	Positive for <i>E. coli</i>

- Look for fluorescence with a 6-watt, 365 nm, UV light within 5 inches of the sample, in a dark environment. Face light away from your eyes and towards the sample.
- If the sample is pink or orange, incubate further (up to 48 hours maximum). A sample that remains pink or orange at 48 hours is negative for coliforms and *E. coli*.
- Samples are negative if at any time after 24 hours there is no red, magenta and/or fluorescence.
- Red/magenta or red/magenta with fluorescence observed before 24 hours is a valid positive. However, after 48 hours, heterotrophs may overwhelm Colisure's inhibition system. Therefore, red/magenta or red/magenta with fluorescence first observed after 48 hours is not a valid positive.

Procedural Notes

- This insert may not reflect your local regulations. For compliance testing, be sure to follow appropriate regulatory procedures. For example, samples run in other countries are incubated at 36±2°C for 24–48 hours.
- Colisure can be run in any multiple tube format. *Standard Methods for the Examination of Water and Wastewater*¹ MPN tables should be used to find Most Probable Numbers (MPNs).
- If sample dilutions are made, multiply the MPN value by the dilution factor to obtain the proper quantitative result.
- Use only sterile, nonbuffered, oxidant-free water for dilutions.
- Colisure is a primary water test. Colisure performance characteristics do not apply to samples altered by any pre-enrichment or concentration.
- Aseptic technique should always be followed when using Colisure. Dispose of in accordance with Good Laboratory Practices.

Quality Control Procedures

1. One of the following quality control procedures is recommended for each lot of Colisure:
 - A. IDEXX-QC Coliform and *E. coli*²: *Escherichia coli*, *Klebsiella variicola*³, and *Pseudomonas aeruginosa*.
 - B. Quanti-Cult³: *Escherichia coli*, *Klebsiella pneumoniae* and *Pseudomonas aeruginosa*.
 - C. Fill three sterile vessels with 100 mL of sterile non-buffered oxidant-free water and inoculate with a sterile loop of ATCC⁴ strains. *Escherichia coli* ATCC 25922/WDCM 00013 or ATCC 11775/WDCM 00090, *Klebsiella variicola*² ATCC 31488/WDCM 00206 and *Pseudomonas aeruginosa* ATCC 10145/WDCM 00024 or ATCC 27853.
2. Follow the P/A Procedure or Quanti-Tray Enumeration Procedure above.
3. Results should match the Result Interpretation table above.

1. Eaton, AD, Clesceri, LS, Greenberg, AE, Rice, EN. *Standard Methods for the Examination of Water and Wastewater* American Public Health Association, 2005. Washington, DC.

2. IDEXX-QC Coliform and *E. coli*—IDEXX Catalog #UN3373-WQC-TCEC

3. Quanti-Cult cultures—IDEXX catalog # WKIT-1001

4. American Type Culture Collection 1-800-638-6597 atcc.org

‡. *Klebsiella pneumoniae* (ATCC 31488/WDCM 00206) has been renamed to *Klebsiella variicola*.

*Colisure, Defined Substrate Technology, DST and Quanti-Tray are trademarks or registered trademarks of IDEXX Laboratories, Inc. or its affiliates in the United States and/or other countries. Quanti-Cult is a trademark or registered trademark of Remel Inc.

Patent information: idexx.com/patents.

© 2019 IDEXX Laboratories, Inc. All rights reserved.

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© 2019 IDEXX Laboratories, Inc. Wszelkie prawa zastrzeżone.